


How to fill a FluoroVette™



Hold the FluoroVette as shown using a 20-200uL pipetter and standard tip

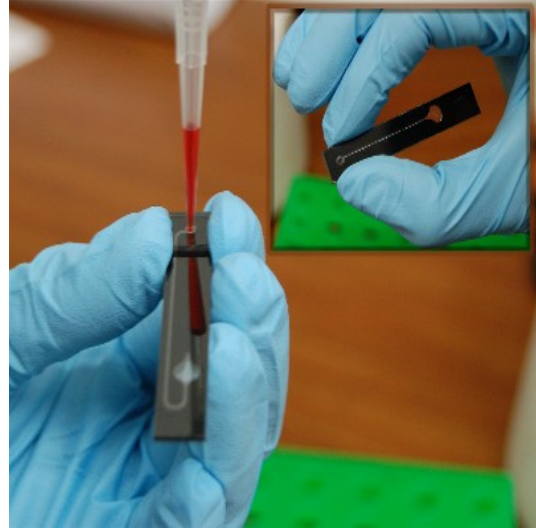

Insert pipette tip at 90° to the pipette interface, Do not obstruct rear outlet



Slowly dispense sample into FluoroVette
The FluoroVette sample volume is 50µL



Bubble trap properly removes air deposits



FluoroVette ready for measurement
With the inlet side of the FluoroVette facing the excitation side of the adaptor



Insert FluoroVette into universal adapter
With the inlet side of the FluoroVette facing the excitation side of the adaptor